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# Thank you,

EN

for purchasing this SCAT Europe product. Please read this operating manual carefully before use and especially observe the safety notices! For service and support, please contact our customer service or your local distributor. For a distributor in your country, please contact us on [www.scat-europe.com](http://www.scat-europe.com).

## Canister

Thank you for purchasing our **electrostatic conductive waste canister**. This waste canister has been designed in such a way to comply with the provisions of CENELEC 50404 (Code of practice for the avoidance of hazards due to static electricity) and the German guideline TRGS 727 (formerly TRBS2153) TRGS 727 can be consulted and downloaded on the website [www.baua.de](http://www.baua.de). Where to obtain CLC/TR50404 in your country, please request at [info@cencenelec.eu](mailto:info@cencenelec.eu)

### **Important notice:**

In order to ensure safe operation we strongly recommend to read this operating manual through carefully before putting your waste canister into service and to comply with the advices when using it. Please check the chemical resistance of the waste canister against the used substances. Also observe the regulations concerning environmental protection and safety in force locally. Please also verify the movability of the floater.

# Content

EN

Overview .....	12
Features.....	13
Proper use.....	14
Grounding .....	15
Safety .....	16
Maintenance and service .....	16

# Overview

EN

# Features

EN

The waste canisters are made of electrostatic conductive plastic (high-density polyethylene PE-HD) and are produced due to quality management standard ISO 9001. The waste canisters exhibit the following electrostatic properties: The waste canisters (inside and outside) are electrostatic conductive.

Fig.	Part No.	Level control	Contents	Thread	UN approval
<b>A</b>	108 421	Viewing strip	10 L	S 50	●
<b>B</b>	108 042	mechanical	10 L	S 60 / 61	-
<b>C</b>	107 953	-	10 L	S 60 / 61	●
<b>D</b>	108 043	mechanical	20 L	S 60 / 61	-
<b>E</b>	108 027	-	20 L	S 60 / 61	●
<b>F</b>	108 193	mechanical	30 L	S 60 / 61	-
<b>G</b>	108 192	-	30 L	S 60 / 61	●
<b>H</b>	107 740	mechanical	60 L	S 70 / 71	-
<b>I</b>	107 940	-	60 L	S 70 / 71	●
<b>J</b>	108 420	mechanical	10 L	S 90	●

**Canisters with UN approval are permitted for transport of hazardous goods on roads and plant premises.**

## Proper use

EN

This electrostatic conductive waste canister is intended for the disposal of chemical substances in the form of liquids or suspensions. The chemicals could be either electrostatic conductive, anti-static or non-conductive. Any use other than the stated agreed use is not permissible.

**Please also note the advice concerning correct grounding.**

**If the product is technically modified by the customer, SCAT Europe GmbH does not assume any liability and guarantee.**

Please pay attention to the chemical resistance of this waste canister to the used chemicals and solvents. According to the wide range of different chemicals we do not assume any liability and guarantee to the compatibility of the waste canister to all chemicals and solvents. For SCAT Europe products the latest most durable materials were selected with special attention to the requirements of working with HPLC systems.

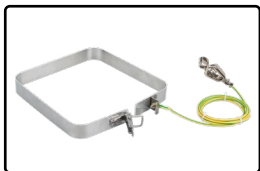
# Grounding

EN

**Important security notice for grounding of the waste canister: In a potentially explosive atmosphere and when working with substances with a low flash-point all objects must be made from electrostatic conductive materials and grounded.**

**The electrostatic conductive waste canisters must therefore be treated like metallic waste canisters and grounded during use.**

Ordering information regarding grounding cables and our extended anti-static equipment you will find at [www.scat-europe.com](http://www.scat-europe.com)

Our products for safety grounding of containers are made for the connection to stainless and electric powered systems. The installation has to be executed by **specialized people only!** Please notice the internal security advises of your company.

**If you have any questions, please contact the person in your laboratory with responsibility for safety matters.**

# Safety

EN

**Before using this waste canister, always check the chemical resistance of the PE-HD against the chemicals and solvents used in your laboratory.**

Ask the supplier of your chemicals about the resistance. You can also find information on the plastic raw material manufacturers' Internet pages and on [www.scats-europe.com](http://www.scats-europe.com). Please follow the instructions given there. Use the waste canister only where chemical resistance is very good, i.e. if no damage or only very slight damage may occur after exposure to the chemical effects for more than 30 days.

**Do not use the waste canister under any circumstances if the resistance is not qualified or if the PE-HD is not resistant. In this case damage to the waste canister is expected and therefore a loss of the waste canister properties will occur.**

Use suitable personal protective equipment to protect yourself against splashing and fumes.

**If you have any questions, please contact the person in your laboratory with responsibility for safety matters.**

## Maintenance and service

**After each use of the waste canister and when carrying out routine maintenance:**

The canister must be cleaned after every usage with water or an appropriate organic solvent to make sure that no insulating coating or diffusion adversely affects the conductivity of the surface and the anti-static properties of the waste canister.

The anti-static properties of a waste canister may be adversely affected by the following factors:

- The liquid or the suspension is sticky and insulating. This may cause a insulating coating to build up on the waste canister.
- The diffusion of a insulating liquid (typically insulating solvents) through the plastic material may have a detrimental effect on its properties over a longer period.

### **Important notice:**

Electrostatic characteristics of the PE-HD material may change over a longer period of time. The resistance to the ground has to be checked at least once a year. The resistance to ground may not be higher than  $10^6$  Ohm at a voltage of 500V.

**If after an annual check of the resistance to ground a worsening of the resistance to ground is detected, the waste canister can no longer considered to be electrostatic conductive and has to be replaced.**

### **Maintenance of the level control floater**

If the floater sticks or is unmovable by sticky or viscous fluids, please clean the level control floater with an appropriate solvent (e.g. water or organic solvent).

**SCAT Europe recommends replacing the canisters routinely, at the latest after 5 years.**

(The UNECE commission issued the ADR (European Agreement concerning the international Carriage of dangerous Goods), which requires an exchange of plastic containers every 5 years from the date of their production (chapter 4.1.1.15). This is conductive to your safety, as the quality of many plastics may change due to the environmental conditions.)

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The latest version of User Manual relating to your product is always available for you to download. Please visit [www.scatt-europe.com](http://www.scatt-europe.com) or scan this QR-Code using the camera of your Smartphone/Tablet.

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